




Configuration Wizard Guide PlasmionNet

Content

1. Welcome	2
2. Getting Started	2
3. Configuration Wizard	3
3.1. Welcome Page.....	4
3.2. License	5
3.3. Device Configuration.....	7
3.4. General Settings & Recording.....	8
3.5. User Management	10
3.6. Summary	11
3.7. Execution	12
3.8. Finish	13
3.9. Cancel / Abort	13
4. Troubleshooting.....	14

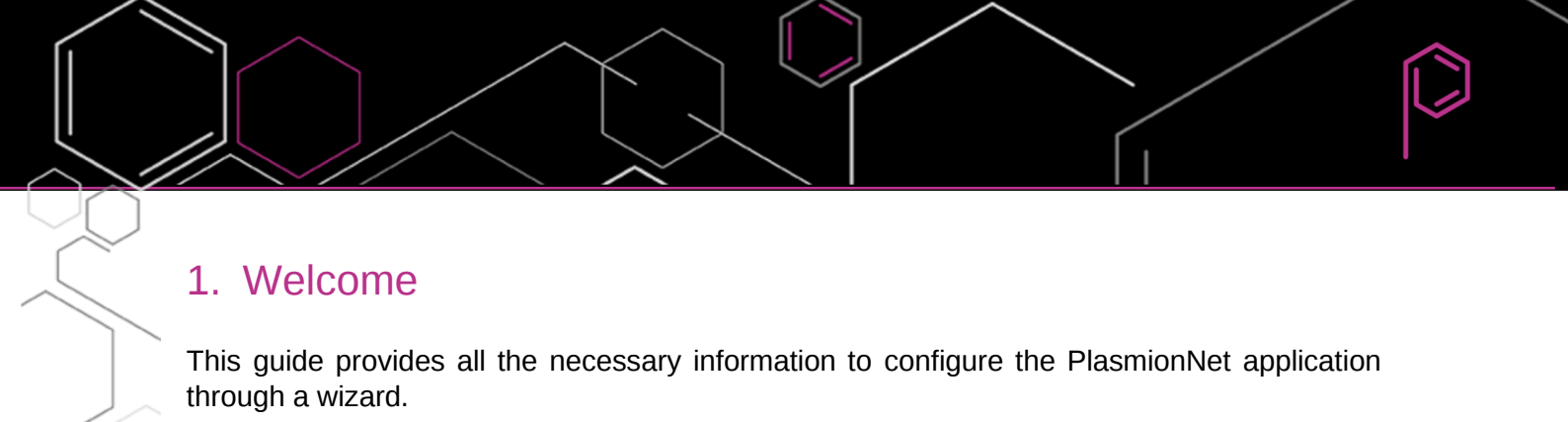

1. Welcome

This guide provides all the necessary information to configure the PlasmionNet application through a wizard.

First, all requirements will be explained. Next, the configuration wizard will be covered. At the end of the document, you will find a section on common errors and troubleshooting.

If you still have any questions or issues, feel free to contact your designated Plasmion representative or use the general contact options.


2. Getting Started

Before starting the configuration process for PlasmionNet, ensure that you meet the following requirements:

- You have successfully installed the PlasmionNet application.
- You have access to the directory where your files shall be stored.

Optional:

If you have already acquired a license file, please have your license file ready for this computer.

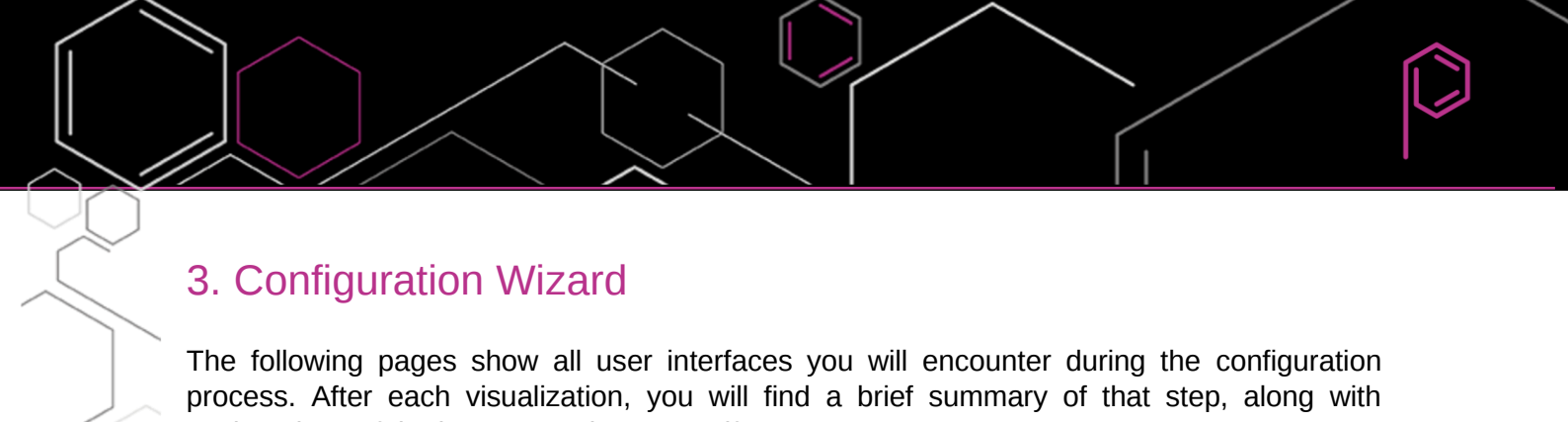


3. Configuration Wizard

The following pages show all user interfaces you will encounter during the configuration process. After each visualization, you will find a brief summary of that step, along with explanations of the buttons on the screen if necessary.

The Wizard will automatically open when you start the Application for the first time on the special file location.

All configuration options from the following *Device Configuration* chapter can also be configured within the application itself. Please refer to the manual for the PlasmionNet application for further instructions.

3.1. Welcome Page

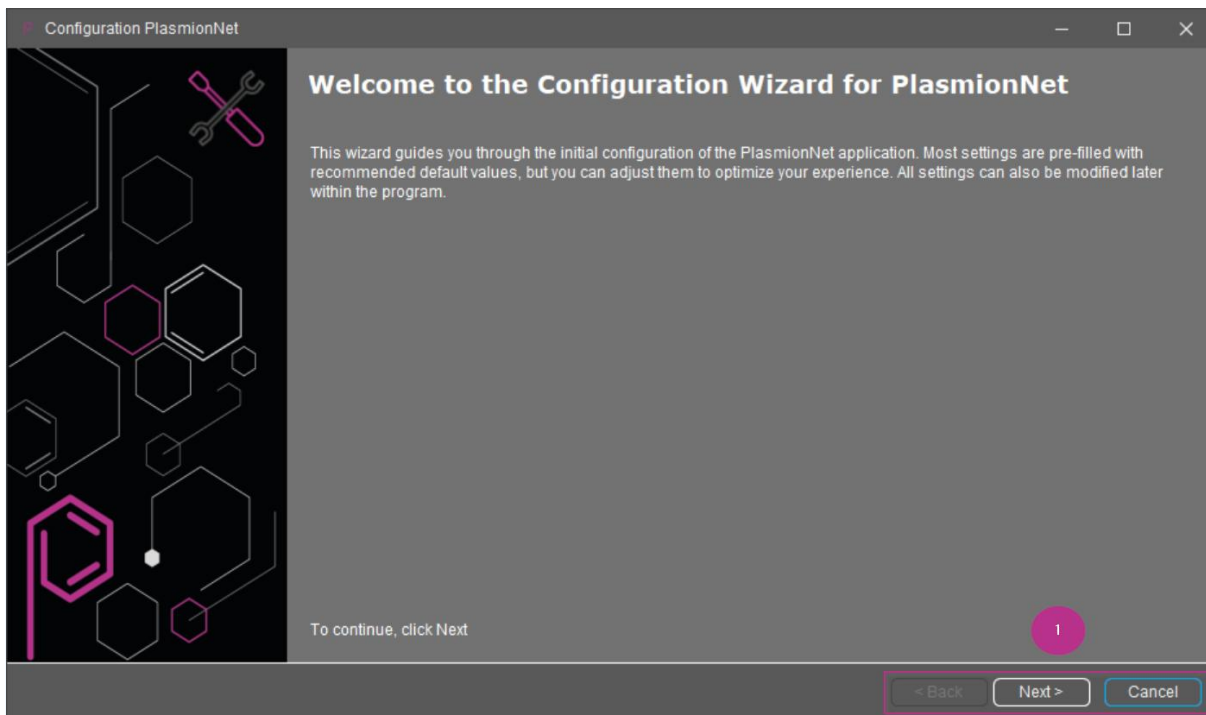

[PR1]

Figure 1 - Welcome Page

The Welcome page provides a brief overview of the actions performed during the execution of this wizard.

In the bottom right corner, you will find the navigation buttons “**Back**”, “**Next**” and “**Cancel**” (marked as **number 1**^[PR2]), which allow you to navigate through the pages of the Wizard.

- The “**Back**” button navigates you to the previous page
- The “**Next**” button validates your inputs and navigates you to the next page if they are valid. Otherwise, a message box will inform you of the validation error.
- The “**Cancel**” button starts the abort process. (For more details, see [Cancel / Abort](#)).

3.2. License

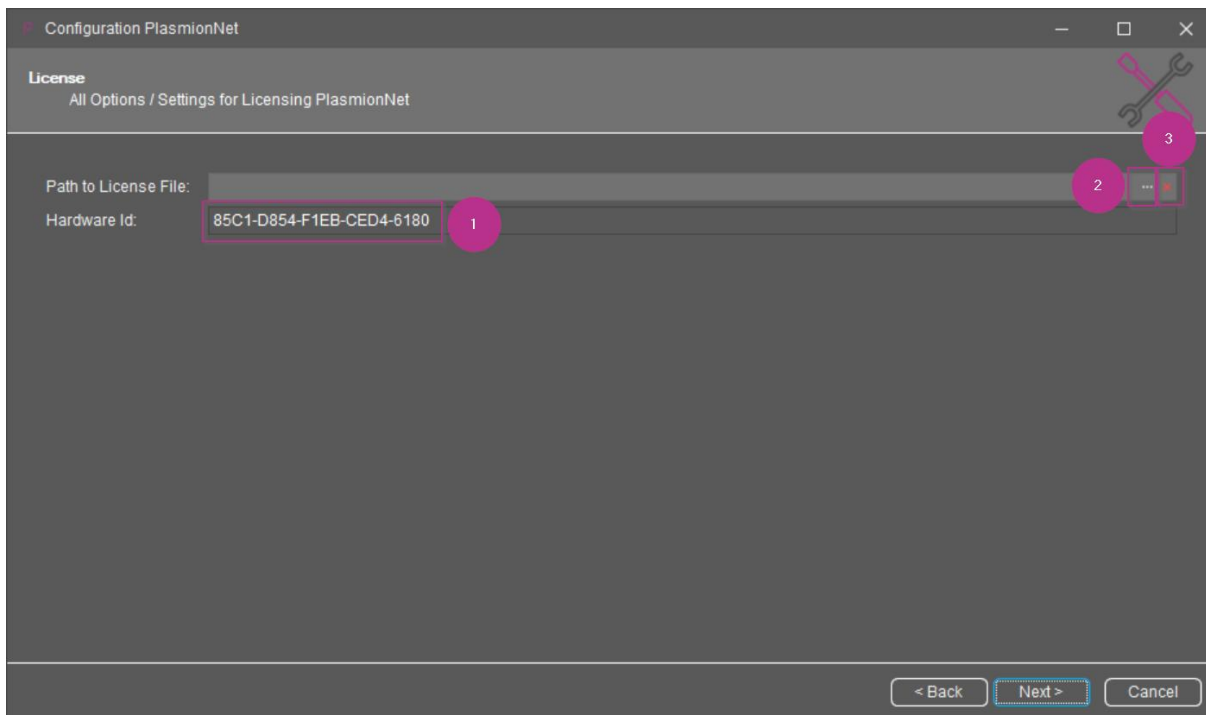


Figure 2 - License & Hardware Id

The License pages collect all information about the licensing of the PlasmionNet Application. You see the Hardware Id (marked as **number 1**) of your computer.

!! Attention: !!

To receive the license file for your computer you must contact your designated Plasmion representative or technical support with the Hardware Id. The representative will then send your license file to you.

With the button (marked as **number 2**) you can now select the license file on your computer (see [Figure 3 - File Selection for License](#)).

The button (marked as **number 3**) allows you to remove the current selection of the file.

!!Attention: !!

The selected license file will be moved from the chosen location to the correct location for the PlasmionNet Application. It is recommended to keep a backup of your license file in case you uninstall the application from your computer.

If you proceed without a license file, a warning will inform you that the application cannot start without a valid license file (see [Figure 4 - Warning if no License is selected](#)).

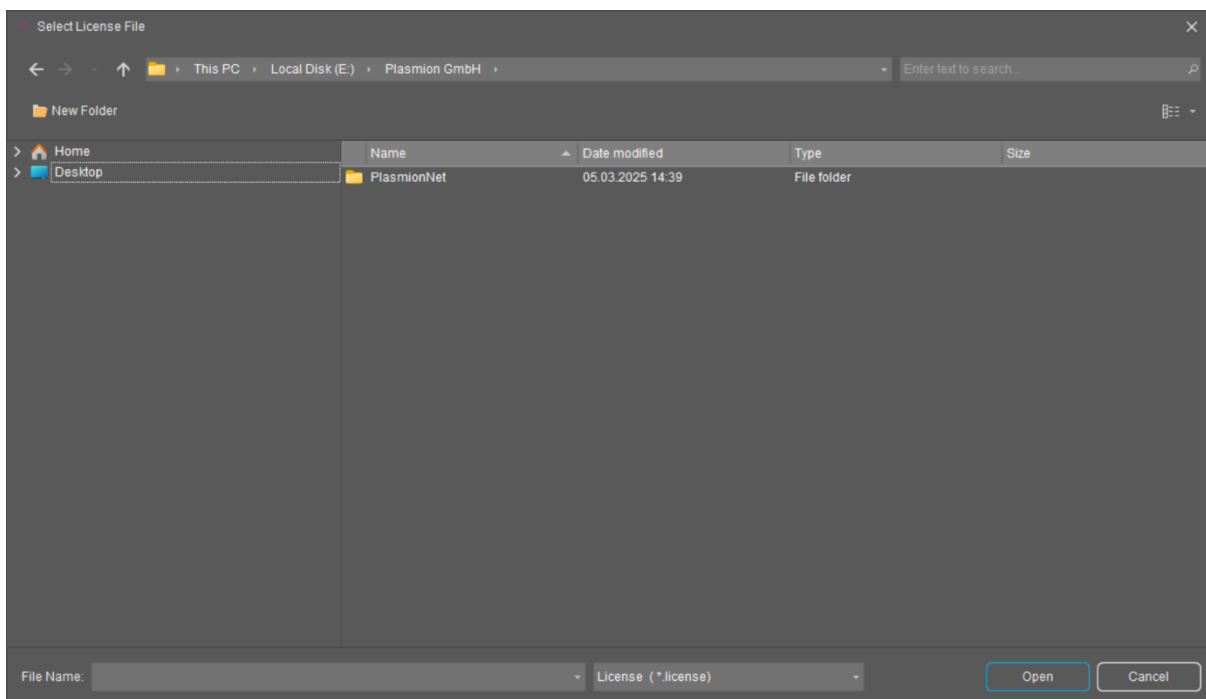


Figure 3 - File Selection for License

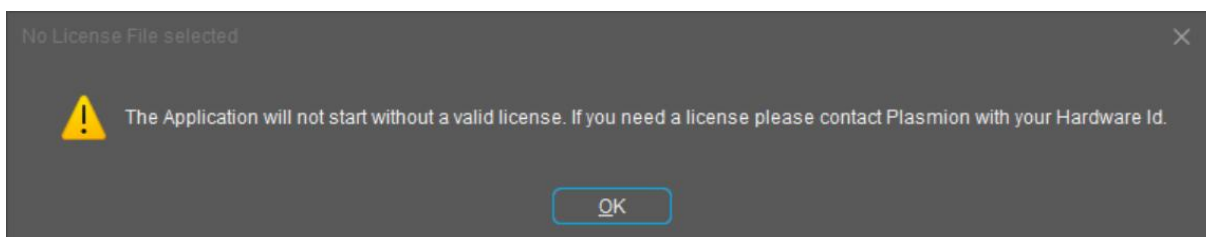


Figure 4 - Warning if no License is selected

3.3. Device Configuration

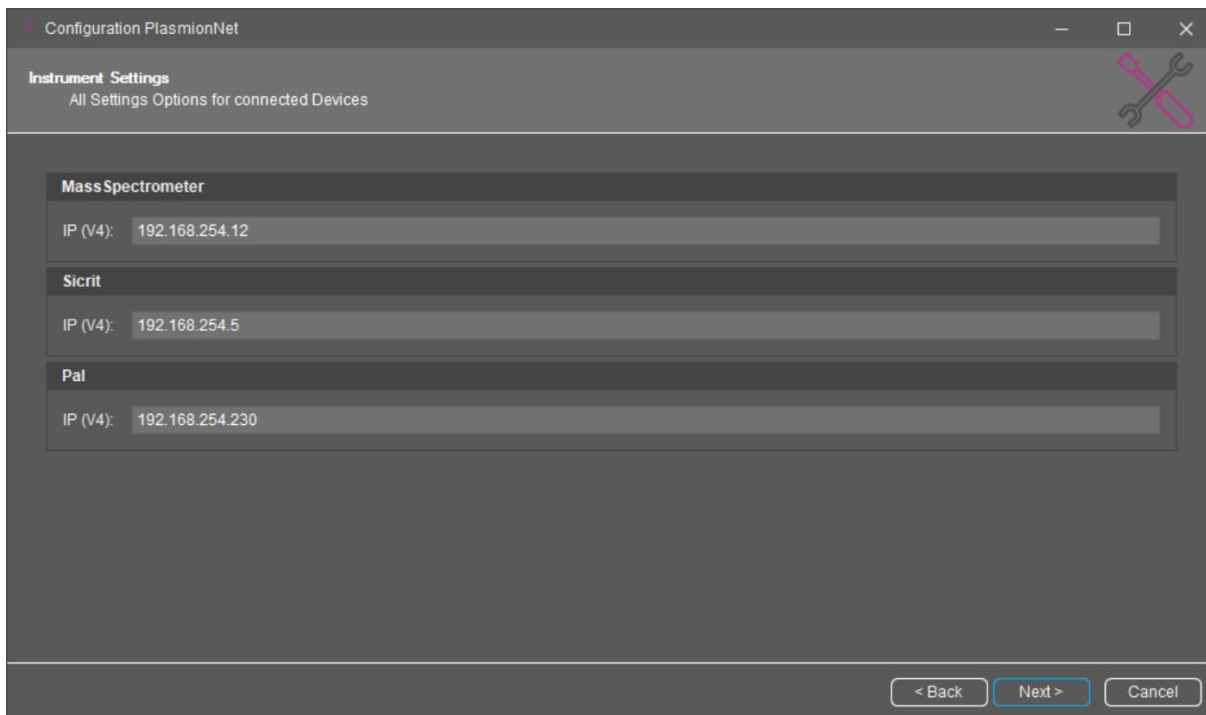


Figure 5 - Connection Parameter for Devices

On this page, you can see the default configuration parameters for the most commonly used devices in the application. You can adjust the settings to match your environment.

When you press the **"Next"** button, the wizard will verify the configured connections to the devices and display a message (see [Figure 6 - Results Connection Test](#)) listing any devices that could not be connected.

Note:

If you have ordered the "HaVoc" product, the values are pre-configured to match the environment.

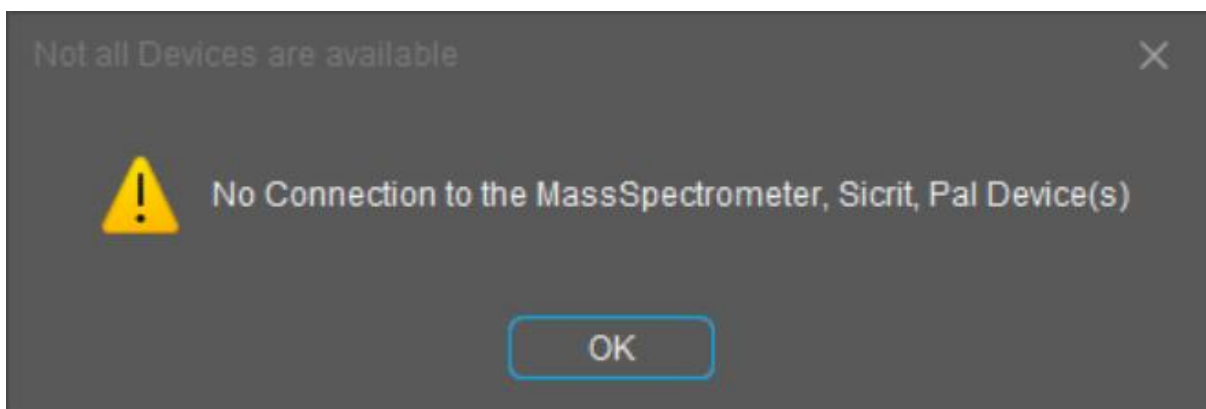


Figure 6 - Results Connection Test

3.4. General Settings & Recording

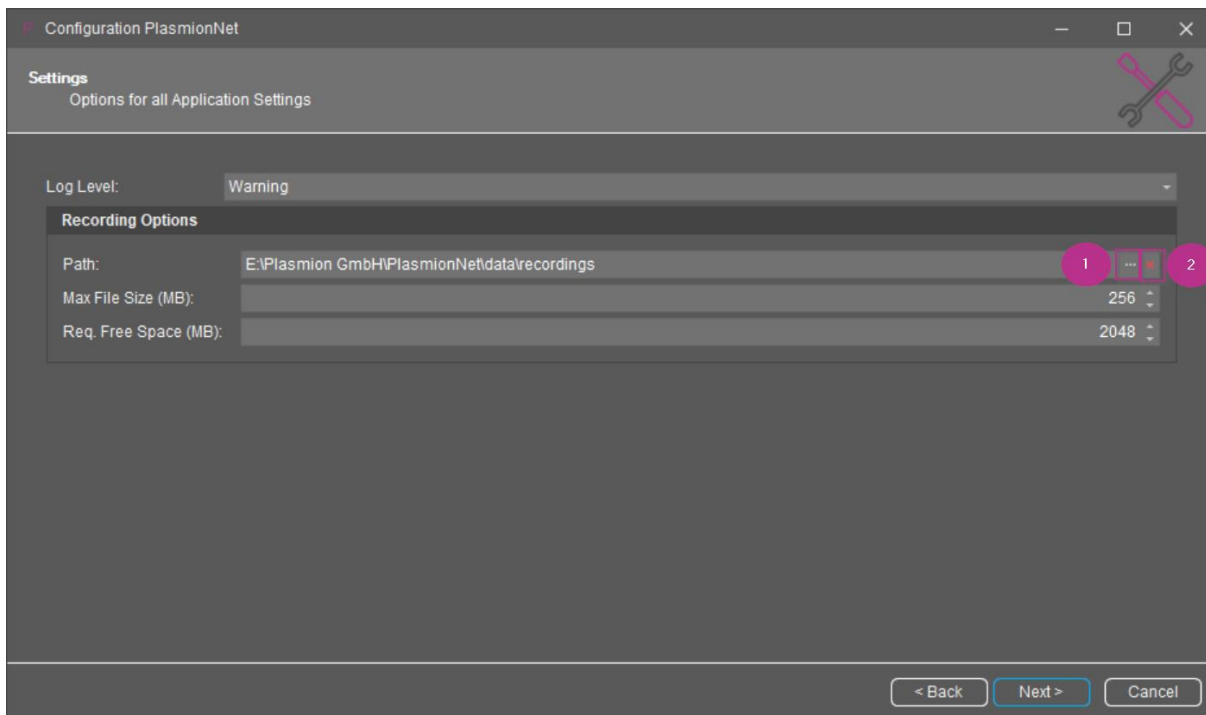


Figure 7 - Recording & General Settings

This page contains all general application settings and a section for recording settings.

The "**Log Level**" setting defines what types of messages are written to the generated log file. Normally, the "**Warning**" level is sufficient. If you encounter frequent issues, you can increase the level to "**Info**" and contact technical support.

The recording settings determine how the application saves measurement results to the file system.

The "**Path**" specifies the directory where all measurement recording files are stored. To select a directory, press the button marked as **number 1**, which will open the selection dialog. To clear the directory, press the button marked as **number 2**. See [Figure 8 - Recording Path Selection](#).

The "**Max File Size (MB)**" setting defines the maximum file size before the recording is split into multiple files. An index marker is added to the file name.

The "**Req. Free Space (MB)**" setting specifies the minimum required free space to ensure recording can proceed. If the available space on the selected drive falls below this limit, you will receive a warning before starting a recording operation in the application.

!!Attention: !!

It is recommended to change the recording path to a separate file location. Otherwise, all recordings will be deleted when you uninstall the application.

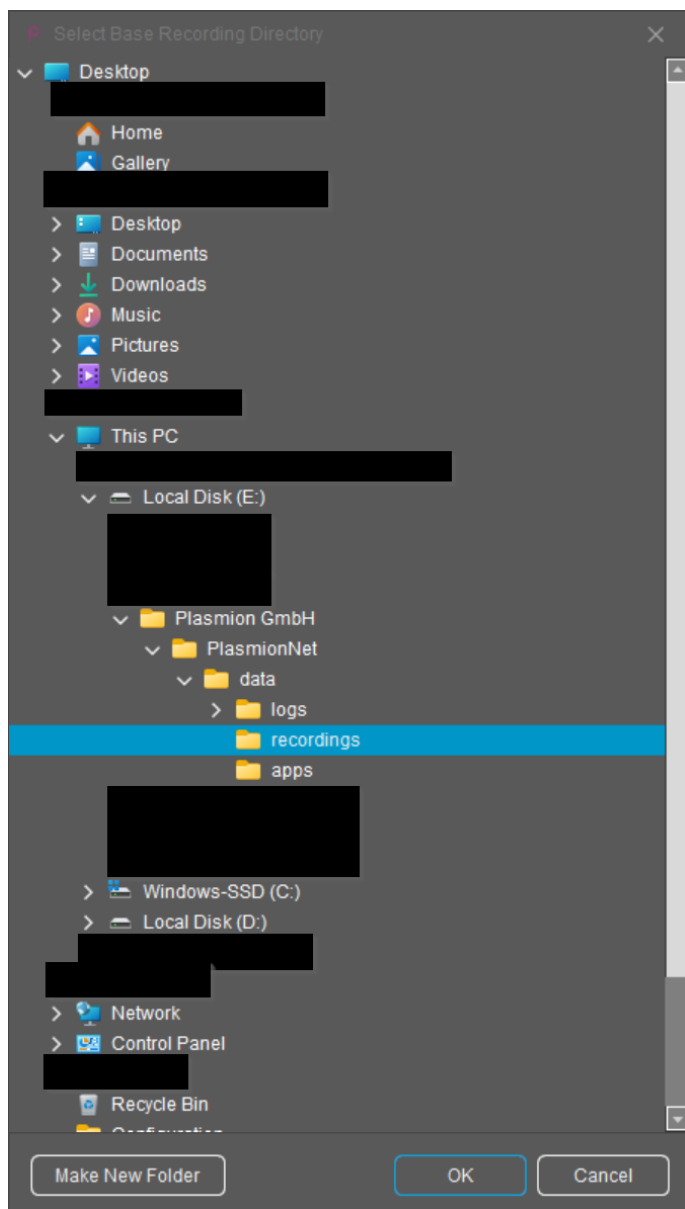


Figure 8 - Recording Path Selection

3.5. User Management

Configuration PlasmionNet

User Management
Configuration for the Application Users

Administrator

Name:

Password:

Password (Confirm):

Operator

Name:

Password:

Password (Confirm):

Remarks: Further Users can create in the "User Management" Area under the Menu Point "Service"

< Back Next > Cancel

Figure 9 - Create default Users

On the User Management page of the wizard, you can configure the default user credentials for the application. You define an **"Administrator"** and an **"Operator"** user.

The **"Administrator"** user has the authority to manage applications, configure the application, and set up users for the application. They are also able to configure settings for a specific application.

The user is allowed to log in to the application and start or execute an application.

The requirements for the new users are:

- A valid name (at least 1 character)
- Both users must have different names
- Password & Password Confirm must match

If any criteria are not met, you will receive a message box with the validation message (see [Figure 10 - Users Requirements Validation](#)).

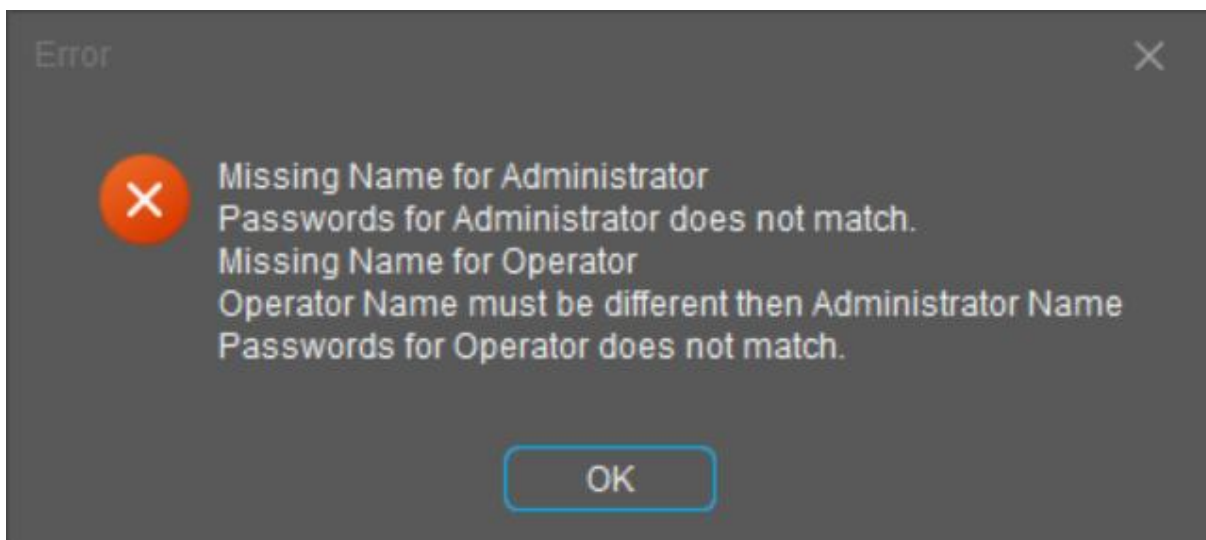


Figure 10 - Users Requirements Validation

3.6. Summary

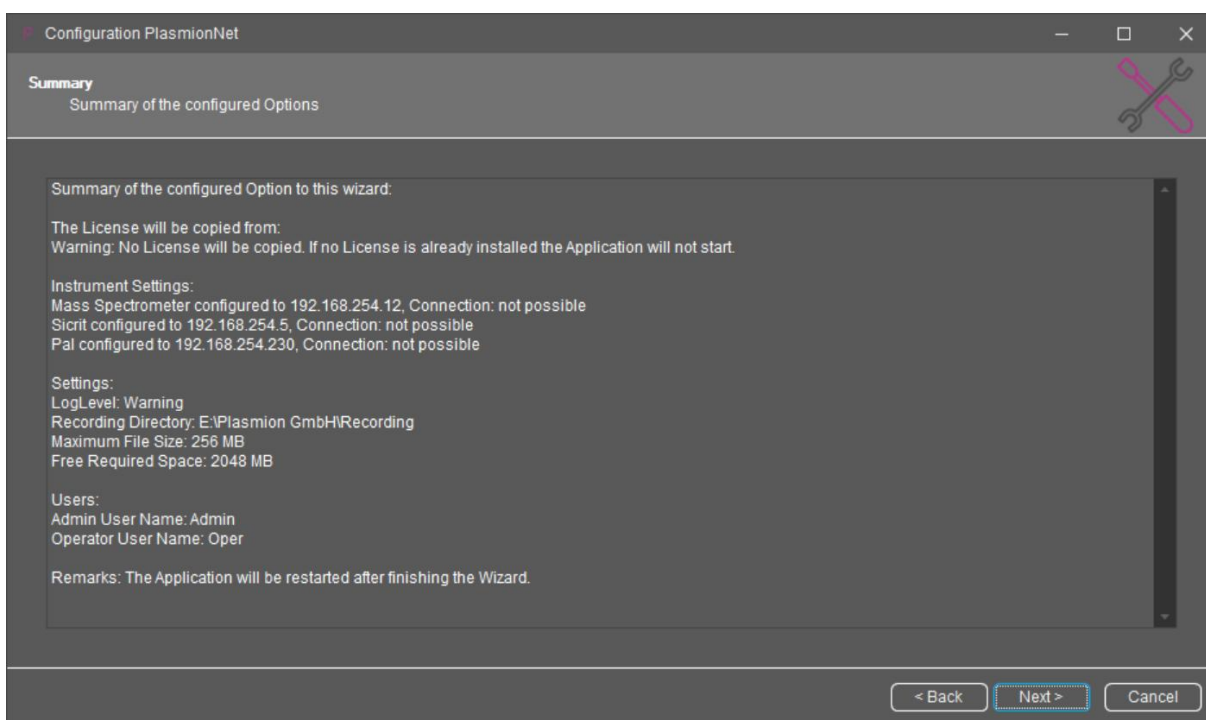


Figure 11 - Generated Summary

The summary page is the final step before the configured inputs are applied. It displays all configured values and provides a summary of the individual pages. It also includes a note that the application will be restarted after successful execution.

3.7. Execution

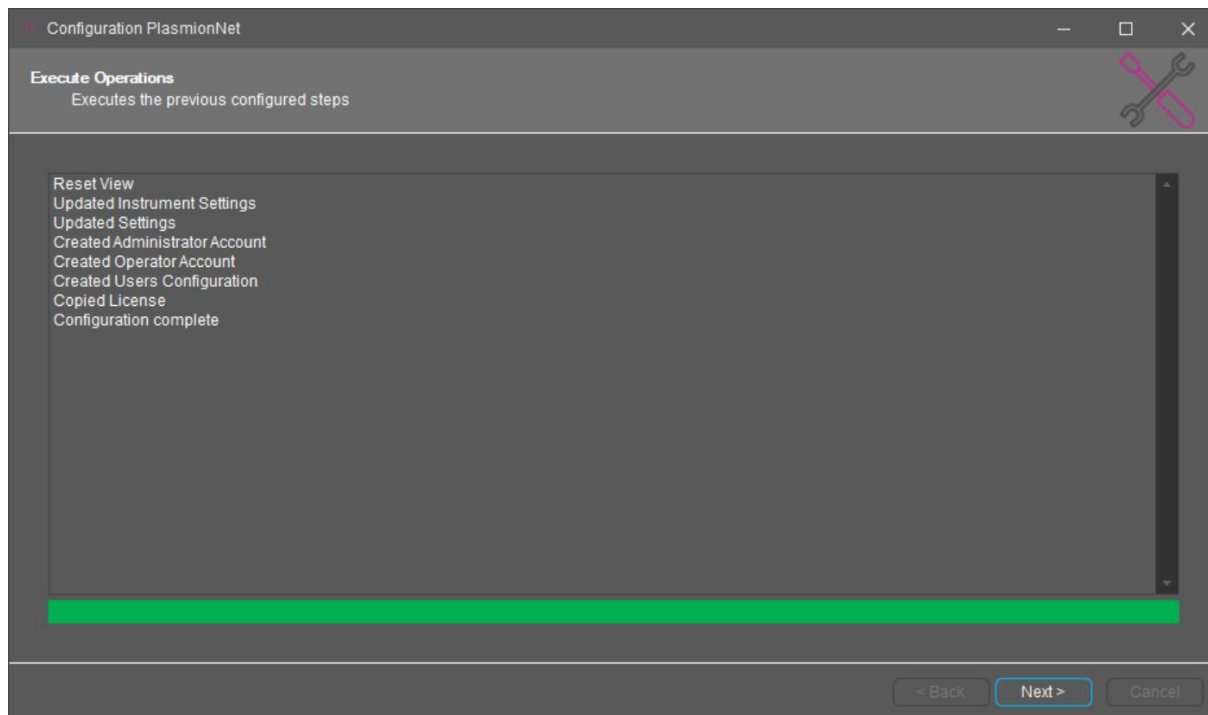


Figure 12 - Wizard execution

This page of wizard shows all the steps being executed. The currently active steps are listed at the end. The progress bar indicates the position of each step in relation to the entire configuration process.

If an error occurs, it will be displayed in the log above the progress bar.

3.8. Finish

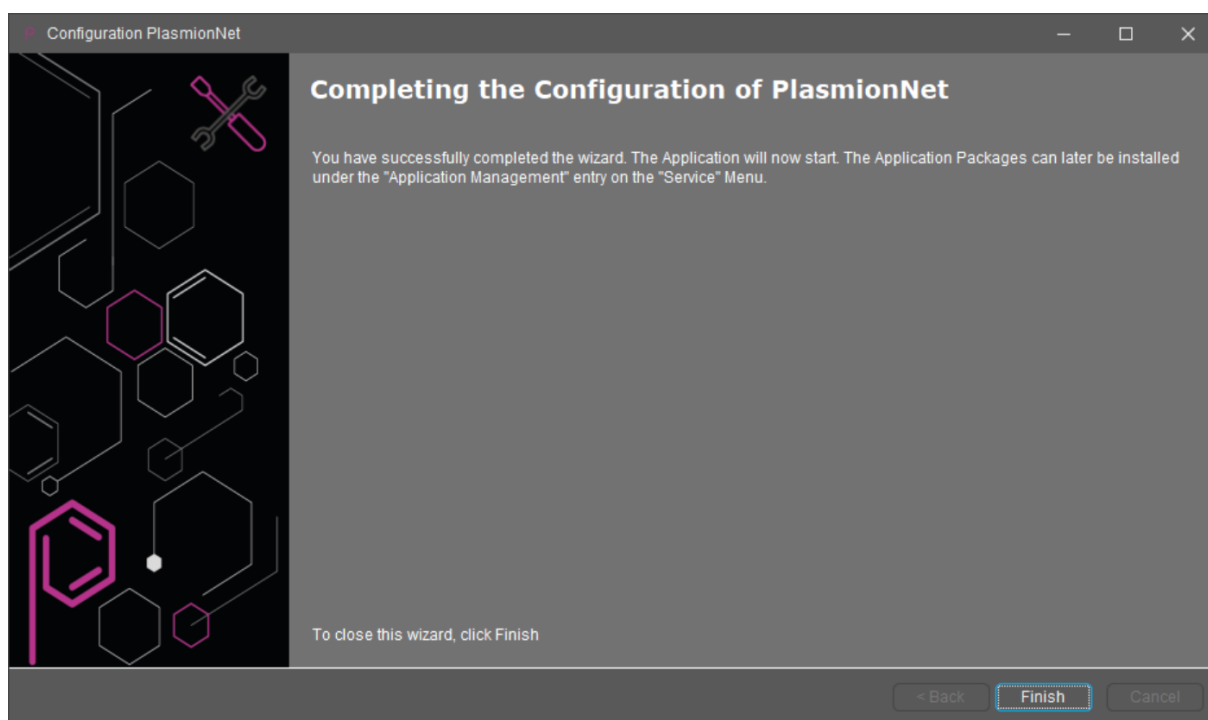


Figure 13 - End of Wizard

This is the final page of the configuration wizard. It indicates that all steps have been successfully configured, and the application will now be restarted with the new values.

3.9. Cancel / Abort

After pressing the "**Cancel**" button on any page of the wizard, you will receive a confirmation message (see [Figure 14 - Cancellation Request](#)). If you confirm the request, the wizard will close without saving any progress. If you deny the request, the wizard will continue on the current page.

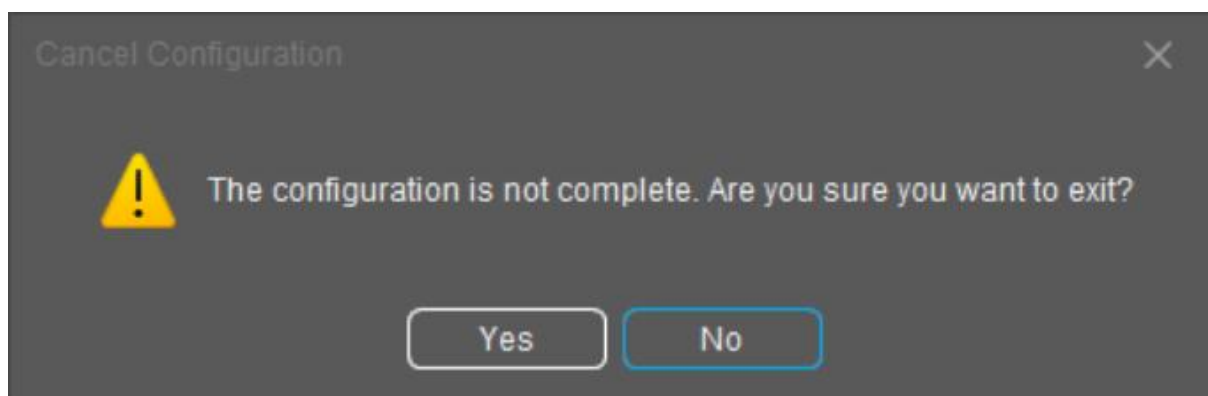


Figure 14 - Cancellation Request

4. Troubleshooting

Problem:

This request is shown when starting the application.

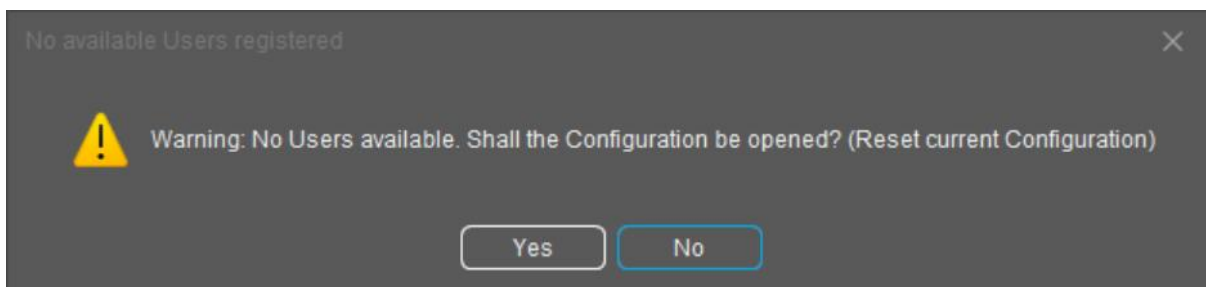


Figure 15 - Request to open the Wizard

Solution:

The application could not find the user configuration file. Please confirm the request to restart the configuration wizard.